

## Slide Staining Rack

### OPERATION

Slide Rack is adjustable to fit tanks up to 22" inside and is adjustable to fit smaller tanks or sinks.

The adjustment is made by sliding the clips to fit the tank either inside or outside.

After several adjustments are made the clips may become loose on the rods. This condition can be corrected by bending the clip with light pressure.

The slide rack is constructed of stainless steel and will resist corrosion under normal use.



Manufactured by:



201 Wolf Drive • PO Box 87 • Thorofare, NJ 08086 • Phone: 1-856-686-1600 • Fax: 856-686-1601

E-mail: [troemner@troemner.com](mailto:troemner@troemner.com) • [www.troemner.com](http://www.troemner.com)

554004-00 (REV 2 - 8/14)